

User instructions

Stereo microscope

KERN

OSE-4

OSE 410, OSE 411, OSE 413
OSE 414, OSE 416, OSE 417

Version 1.0
01/2015





KERN OSE-4

Version 1.0 01/2015

User instructions

Stereo microscope

Table of contents

1	Before use	3
1.1	General notes.....	3
1.2	Notes on the electrical system.....	3
1.3	Storage	4
1.4	Maintenance and cleaning.....	5
2	Nomenclature	6
3	Basic data.....	7
4	Assembly	8
5	Operation and functionality.....	10
5.1	Getting started	10
5.2	Adjusting the interpupillary distance	10
5.3	Adjusting the magnification	10
5.4	Dioptre adjustment and focussing	11
5.5	Adjusting the stand	12
5.6	Using eye cups / High Eye Point eyepieces	13
5.7	Lighting control	14
5.8	Using external illumination units.....	15
5.9	Changing the bulb	15
6	Optical data	16
7	Features.....	16
8	Trouble shooting.....	17
9	Service	18
10	Disposal.....	18
11	Further information.....	18

1 Before use

1.1 General notes

You must open the packaging carefully, to make sure that none of the accessories in the packaging fall on the floor and get broken.

In general, microscopes should always be handled carefully because they are sensitive precision instruments. When using or transporting the microscope it is particularly important to avoid abrupt movements, as this may damage the optical components.

You should also avoid getting dirt or finger prints on the lens surface, because in most cases this will reduce image clarity.

To maintain the performance of the microscope, it must never be disassembled. So components such as lenses and other optical elements should be left as they were before use. Also the electrical parts in the base of the device must not be tampered with, as in this area there is an additional risk of triggering an electric shock.

1.2 Notes on the electrical system

(OSE 413, OSE 414, OSE 416, OSE 417)

Before connecting to a mains power supply, you must make sure that you are using the correct input voltage. The mains cable supplied has an external power supply. The necessary values are shown on this power supply. If you do not comply with these specifications, then fires or other damage to the device could occur.

The main switch must also be switched off before the mains cable is connected. In this way you will avoid triggering an electric shock.

If you are using an extension cable, then the mains cable you use must be earthed.

When carrying out any procedures whereby you come into contact with the electrical system of the device, such as, for example, changing the bulb, only carry out these procedures when the power is disconnected.

1.3 Storage

You should ensure that the device is not exposed to direct sunlight, temperatures which are too high or too low, vibrations, dust or a high level of humidity.

The ideal temperature range is between 0 and 40°C and a relative humidity of 85% should not be exceeded.

The device should always be located on a rigid, smooth, horizontal surface.

For devices with pillar stands, the microscope holder must not be rotated back too far. If you do this, there is a risk that the microscope could tip over.

When the microscope is not being used, you should fit the objective cap and cover the microscope with the enclosed dust protective cover.

If the eyepieces are being stored separately, the protective caps must be fitted to the tube connectors. In most cases, if dust and dirt gets inside the optical unit of a microscope this can cause irreversible errors or damage.

The best way to store accessories which consist of optical elements, such as, for example, eyepieces and objectives, is in a dry box with desiccant.

1.4 Maintenance and cleaning

In any event, the device must be kept clean and dusted regularly.

If any moisture should occur, before you wipe down the device you must ensure that the mains power is switched off.

When glass components become dirty, the best way to clean them is to wipe them gently with a lint-free cloth.

To wipe oil stains or finger prints off the lens surface, moisten the lint free cloth with a mixture of ether and alcohol (70 / 30 ratio) and use this to clean the lens.

You must be careful when handling ether and alcohol, as these are highly flammable substances. You must therefore keep it away from naked flames and electrical devices which can be switched on and off, and only use it in well-ventilated rooms.

However organic solutions of this type should not be used to clean other components of the device. This could lead to damage to the paint finish. To do this, it is sufficient to use a neutral cleaning product.

You could also use the following cleaning products to clean the optical components:


- Special cleaner for optical lenses
- Special optical cleaning cloths
- Bellows
- Brush

When handled correctly and checked regularly, the microscope should give many years of efficient service.

Should repairs still be necessary, please contact your KERN dealer or our Technical Department.

2 Nomenclature

Using the OSE 416 as an example
(with reflected and transmitted light illumination)

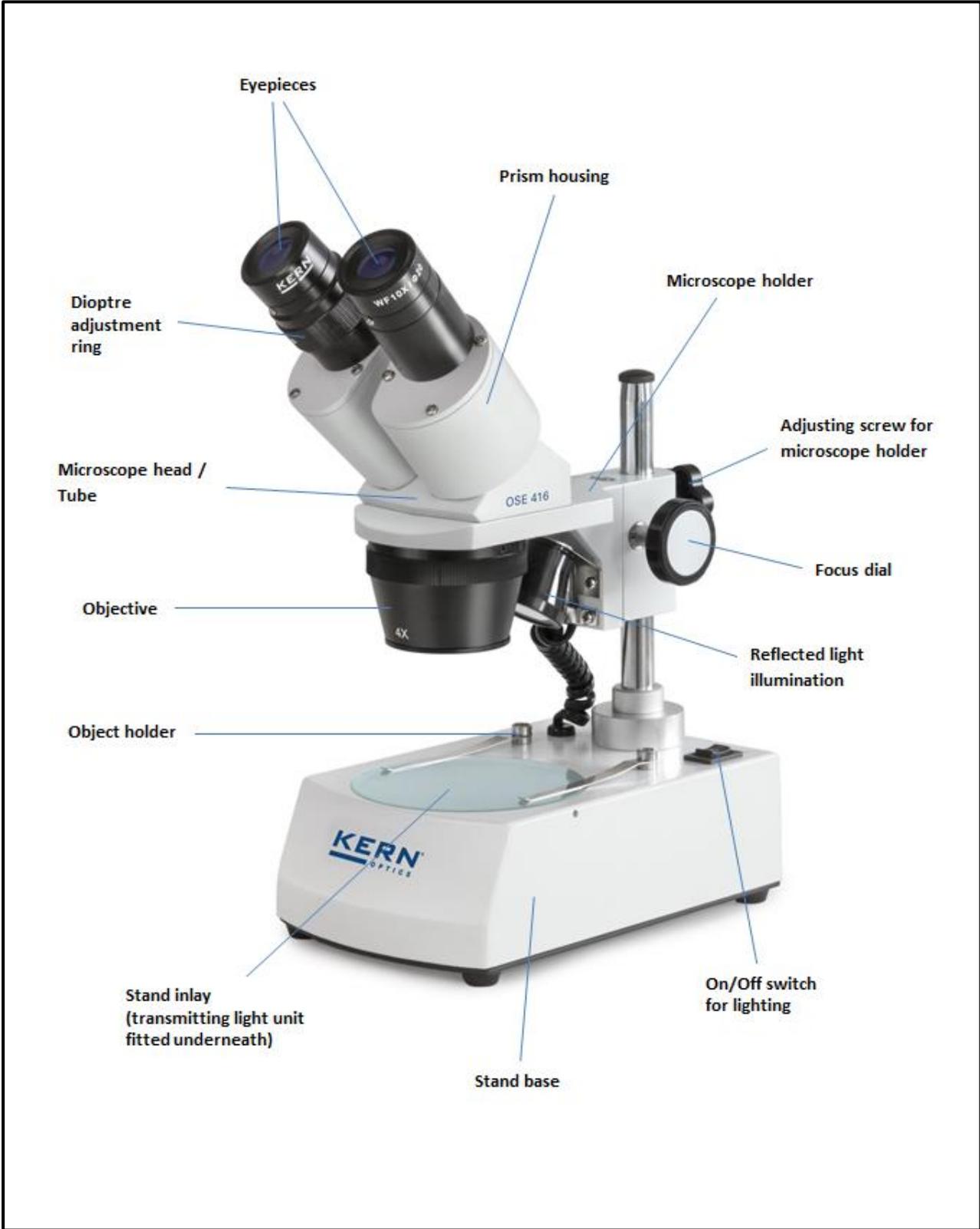

3 Basic data

Optical system	Greenough
Dimmable lighting	No
Tube	angled at 45°
Interpupillary distance	55 – 75 mm
Dioptre adjustment	One-sided (left)
Packaging dimensions WxDxH	195 x 162 x 345 mm

Standard configuration

Model	Tube	Eyepiece	Field of view	Objective	Stand	Illumination
KERN			mm			
OSE 410	Binocular	WF 10x Ø 20 mm	Ø 20	1x / 3x	Pillar	-
OSE 411	Binocular	WF 10x Ø 20 mm	Ø 20	2x / 4x	Pillar	-
OSE 413	Binocular	WF 10x Ø 20 mm	Ø 20	1x / 3x	Pillar	0.21W LED (reflected light)
OSE 414	Binocular	WF 10x Ø 20 mm	Ø 20	2x / 4x	Pillar	0.21W LED (reflected light)
OSE 416	Binocular	WF 10x Ø 20 mm	Ø 20	1x / 3x	Pillar	0.21W LED (reflected light); 0.21W LED (transmitted light)
OSE 417	Binocular	WF 10x Ø 20 mm	Ø 20	2x / 4x	Pillar	0.21W LED (reflected light); 0.21W LED (transmitted light)

4 Assembly

The first step is to position the **microscope stand on a firm, level surface**.

The holder with the head is already on pillar of the stand, but you must always check whether it is **fixed** securely and that it is in the **correct position** (preferably centrally and facing to the front).

Please see section 5.5 for more details on adjusting the stand.

Ideally, the holder and head are then parallel on the central axis of the stand base (see figure on page 9).

Both eyepieces are already fitted in the tube. You just need to remove the protective film which covers them.

When doing this make sure that you **do not touch the optical lenses with your fingers**.

In general you should avoid using two eyepieces with different magnifications.

When **changing the eyepieces**, you must also loosen the **small silver screw** under the eyepiece on the tube connectors and once the new eyepiece is fitted, then it must be re-tightened.

With regard to the use of a device **with transmitted light illumination** (OSE 416, OSE 417) you must make sure that the **frosted glass stand insert is fitted in the centre of the stand base**, so that the transmitted light can be used correctly.

Additional optional attachments:

- The eye cups supplied with the microscopes can be fitted to the eyepieces. (see section 5.6)

5.4 Dioptre adjustment and focussing

A special feature of stereo microscopes is that they are fitted with an optical unit which has a relatively high depth of field. In order to be able get the most benefit from this feature, each user must synchronise the focussing mechanisms for themselves.

The steps to do this are described in the following section.

1. Place the object to be observed on the surface under the objective.
2. Use the objective to set the smallest possible magnification.
3. Look through the righthand eyepiece (without dioptre adjustment ring) with the right eye and bring the object into focus by using the focus control dials.
4. Now set the largest magnification.
5. Once again, still only looking through the righthand eyepiece, bring the object into focus
6. Then set the smallest possible zoom factor again.
7. Then look through the lefthand eyepiece with the left eye (with dioptre compensation ring) and bring the object into focus, by turning the dioptre compensation ring left or right at the right point.
8. In order to get the highest level of accuracy when adjusting the focus, you should repeat steps 4-7.
9. In this way, the object being observed will be in focus on any magnification setting.

5.5 Adjusting the stand

Torque of the focus wheels

You can adjust the torque of the focus wheels by turning the ring which is fitted along the axis of the lefthand focus wheel.

To do this you need a special wrench which is included with delivery. The ring has holes which the wrench can hook into, so that you can then turn it in the desired direction.

Depending on the direction of the turn, the torque will be increased or decreased.

On one hand, this function can help to make it easier to adjust the focus and on the other hand it can prevent the microscope head from slipping down unintentionally. In this way you can avoid possible damage which could occur if the objective lens and the object being observed should collide.

Height adjustment



Fixing the microscope holder

With a stereo microscope on a pillar stand, the microscope head can be height adjusted using the focus wheels and in addition, the microscope holder can be fixed at any point on the pillar, depending on the application requirements.

To fix the position, use the adjusting screw on the holder (*see figure*).

5.6 Using eye cups / High Eye Point eyepieces

The eye cups supplied with the microscope can basically be used at all times, as they screen out intrusive light, which is reflected from light sources from the environment onto the eyepiece, and the result is better image quality.

But primarily, if eyepieces with a high eye point (particularly suitable for those who wear glasses) are used, then it may also be useful for users who don't wear glasses, to fit the eye cups to the eyepieces.

These special eyepieces are also called High Eye Point eyepieces. They can be identified by the glasses symbol on the side. They are also marked in the item description by an additional "H" (example: HSWF 10x Ø 23 mm).

When fitting the eye cups, make sure that the dioptre setting is not moved. We would therefore advise that you hold the dioptre adjustment ring on an eyepiece with one hand while you fit the eye cup with the other.

Before using the microscope, users who wear glasses must remove the eye cups, which you may find on High Eye Point eyepieces.

As the eye cups are made of rubber, you must be aware that when you are using them, they can become slightly dirty through grease residues. In order to maintain hygiene, we would therefore recommend that you clean the eye cups regularly (e.g. with a damp cloth).



Eye cups



High Eye Point eyepiece
(identified by the glasses symbol)

5.7 Lighting control

(OSE 413, OSE 414, OSE 416, OSE 417)

OSE 413, OSE 414 (reflected light)

To operate the reflected light, there is a main switch on the right, rear corner of the stand base surface which you can use to switch the lighting on and off.

The light intensity cannot be adjusted.

OSE 416, OSE 417 (reflected and transmitted light)

To operate the reflected and transmitted light, there is a main switch on the right, rear corner of the stand base surface.

This switch has three positions:

O position	Lighting is switched off
I position	Transmitted light is switched on
II position	Reflected light is switched on

Reflected light and transmitted light cannot be used at the same time.

The light intensity cannot be adjusted.

5.8 Using external illumination units

If, when using a microscope in its standard version, the lighting is not suitable for the application, then it often makes sense to fit an external lighting unit to overcome this problem.

The lighting units which are suitable for devices of the OSE-4 series, are goose neck lighting units (*see figure*). These are available as LED as well as halogen versions and also have an on/off switch or different controller.



Typical goose neck lighting unit

Using goose neck lighting units

Depending on the requirements, a goose neck lighting unit is placed next to, in front of or behind the microscope. The light source is in the housing of the unit and is emitted through one or several optical fibre cables.

These cables are flexible and therefore offer a large number of positioning options, so that the object being observed is perfectly illuminated.

5.9 Changing the bulb

LED

The devices in the OSE-4 series with lighting are all fitted with LED bulbs.

Due to the long service life of an LED lighting system, for these microscopes it will not be necessary to simply change a bulb.

Problems with the lighting unit would therefore, in most cases, be caused by defects in the electrical system. If this is the case, then our Technical Service will be able to help.

6 Optical data

Eyepiece	Specifications - Objectives				
	Magnification	1x	2x	3x	4x
WF 5x	Total magnification	5x	10x	15x	20x
	Field of view mm	Ø 20	Ø 10	Ø 6,7	Ø 5
WF 10x	Total magnification	10x	20x	30x	40x
	Field of view mm	Ø 20	Ø 10	Ø 6,7	Ø 5
WF 15x	Total magnification	15x	30x	45x	60x
	Field of view mm	Ø 15	Ø 7,5	Ø 5	Ø 3,7
WF 20x	Total magnification	20x	40x	60x	80x
	Field of view mm	Ø 10	Ø 6,5	Ø 4,3	Ø 3,2
Working distance		57 mm	57 mm	57 mm	57 mm

7 Features

Model outfit		Kern model						Order number
		OSE 410	OSE 411	OSE 413	OSE 414	OSE 416	OSE 417	
Eyepieces	WF 5x / Ø 16.2 mm	oo	oo	oo	oo	oo	oo	OZB-A4101
	WF 10x / Ø 20 mm	●●	●●	●●	●●	●●	●●	OZB-A4102
	WF 15x / Ø 15 mm	oo	oo	oo	oo	oo	oo	OZB-A4103
	WF 20x / Ø 10 mm	oo	oo	oo	oo	oo	oo	OZB-A4104
Stand	Pillar, without lighting	●	●					
	Pillar, with 0.21W LED lighting (reflected light)			●	●			
	Pillar, with 0.21W LED lighting (Transmitted light and reflected light)					●	●	
Stand inlay	black-white / Ø59.5 mm	●	●	●	●			OZB-A4816
	Frosted glass / Ø95 mm					●	●	OZB-A4805
	black-white / Ø95 mm					●	●	OZB-A4806
External illumination	For information on external lighting units, please see the Kern Optics main catalogue, from page 68 and visit our website www.kern-sohn.com							

● = Standard configuration

o = Option

8 Trouble shooting

Electrical system

Problem	Possible causes
The lighting unit (if fitted) cannot be switched on	The power cable is either not connected or not connected correctly
	The bulb is not fitted
	The bulb has blown
	The fuse has blown
The bulb has blown	The brightness control is set to the lowest level
	The wrong bulb has been used
The bulb flickers	The input voltage was too high
	The bulb is not correctly fitted
The bulb brightness is not sufficient	The lamp is worn out
	The wrong bulb has been used
	The input voltage is too low

Optical unit

Problem	Possible causes
You can see two images	The gap between the eyes is not set correctly
	The magnifications of the eyepieces do not match
There is dirt in the visual field	There is dirt on the object being observed
	There is dirt on the eyepiece surface
The image is unclear	There is dirt on the objective surface
The focus wheels are jammed	The torque of the focus wheels is set too high
The microscope head slips down while you are viewing the object	The torque of the focus wheels is set too low
Eyes get tired easily	The dioptre adjustment is not correct
	The brightness adjustment is not correct

9 Service

If, after studying the user manual, you still have questions about commissioning or using the microscope, or if unforeseen problems should arise, please get in touch with your dealer. The device may only be opened by trained service engineers who have been authorised by KERN.

10 Disposal

The packaging is made of environmentally-friendly materials, which you can dispose of at your local recycling centre. Disposal of the storage box and device must be carried out by the operator in accordance with all national or regional laws in force in the location of use.

11 Further information

The illustrations may differ slightly from the product.

The descriptions and illustrations in this user manual are **subject to change without notice**. Further developments on the device may lead to these changes.



All language versions contain a non-binding translation.
The original German document is the binding version.